What is biowaste?

Society has long used biowaste in the form of compost for improving or fertilising soils, or for releasing heat by burning. But in the last 20 years or so, there has been growing realisation that biowaste can also be used to generate useful chemicals, by subjecting it to high temperatures in the absence of air (known as pyrolysis)

The term 'biowaste' covers a wide range of materials – from crop stalks to woodchip, and from spent hops to seaweed. But what it all has in common is that it was recently living material that was sequestering carbon from the atmosphere, and so inherently avoids the large CO2 footprint of fossil fuels.

These products can be very varied, says Dr Daniel Nowakowski, an expert in pyrolysis at Aston University, UK: "By heating up biowaste in a carefully controlled fashion, we can generate solids, liquids and gases with a wide range of interesting uses".

Deciding what to do with a feedstock

As Lecturer in Chemical Engineering and Applied Chemistry at Aston's Energy and Bioproducts Research Institute (EBRI), Dr Nowakowski has first-hand knowledge about the types of biowaste sources and the ways in which the resulting products can be used – from carbon-storing soil conditioners to sustainable fuels.

This knowledge is put to use every day, not just through the team's fundamental research, but through their project work for government agencies, organisations and companies looking to do more with their biowaste. When they're given a biowaste sample to investigate, he says, we ask one question: "How can we tailor the pyrolysis process to derive products with the greatest value or benefit?".

For example, he says, they might want to maximise yields of a particular product or eliminate an unwanted contaminant. "Working out the processing conditions that will do this requires experience and research, but also a lot of experimentation: a single Ph.D. student can work for three years on optimising the yield of a single component from a pyrolysed feedstock!".

Pyrolysis – Process and products

Before going further, it's worth first understanding what pyrolysis is. Otherwise known as 'heat treatment', pyrolysis involves raising the temperature of a material in a reactor to 400–600 °C in the absence of oxygen. The material needs to be fairly dry, with typically less than 20% water, to avoid compromising energy efficiency and causing phase separation of pyrolysis liquids. This closely mirrors the age-old process of making charcoal, says Dr Nowakowski, but with much greater control over the conditions. Also, he says, nothing goes to waste: "Unlike in conventional charcoal production, we're not just interested in the solid matter left behind – we also capture all the vapours and gases emitted, in order to extract value from them too", he explains.

During the pyrolysis process, there is extensive rearrangement and degradation of the bonds in the lignin, cellulose and hemicellulose and other compounds that make up a typical biowaste feedstock. This leads to the production of three classes of materials:

- **Biochar** is a highly porous form of carbon, and typically takes the form of a black powder. It's primarily of interest as a soil conditioner that can store carbon in the soil for extended periods, in a process known as carbon sequestration and storage (CCS). There are numerous standards for biochar, and biochar classes have also been recognised on the basis of the carbon content, particle size, and presence of plant nutrients such as potassium and nitrogen. Biochar can also be 'engineered' further into catalyst supports, polymer fillers, and absorbents for heavy metals.
- **Bioliquids** contain a complex mix of carboxylic acids, phenols (and its derivatives), esters, ketones, anhydrides, ethers, and

oxygen-containing heterocycles. During pyrolysis, they're released as vapours, but are subsequently condensed into liquids. The water-miscible (aqueous) fractions, because of their acid content, are also known as 'wood vinegar', which has attracted interest as a fungicide and plant growth enhancer. The water-immiscible ('organic') fractions are rich in aromatic compounds, which can either be separated from each other and used as chemical 'building blocks', or catalytically upgraded to remove the oxygen atoms and give hydrocarbon fuels.

• Pyrolysis gas (also called 'pyrogas') includes methane, ethane and other hydrocarbons up to C5, hydrogen, carbon monoxide and traces of carbon dioxide. Gaseous hydrocarbons can be used to fuel the pyrolysis reactor, while hydrogen and carbon

monoxide (syngas) can be used to produce ammonia or methanol. Hydrogen is also of interest as a 'green' fuel for transport, particularly for heavy goods vehicles (for which electrification is not currently practical).

Tailoring the pyrolysis process

The amounts of biochar, bioliquids and biogases produced vary greatly depending upon the volatile content of the original material and the processing conditions, says Dr Sarah Asplin, a Research Associate at EBRI who is using her knowledge and skills from her Ph.D. in the field of biomass pyrolysis. She explains that slow, relatively low-temperature pyrolysis of biomass in a batch reactor might yield up to 50% biochar and 25% each of bioliquid and biogas. In contrast, quick, hightemperature pyrolysis of high-quality woodchip in a fluidisedbed reactor could result in up to 75% bioliquid, with most of the remainder being gases, and very little biochar.

Along with feedstock quality, all these factors feed in to decisions about what to do with a particular feedstock, says Dr Asplin: "For example, clean feedstocks, especially those that are lignin-rich,

LOW CARBON, BIG IMPACT: MAKING THE MOST OF BIOWASTE

Making useful materials from biological waste streams using pyrolysis

Pyrolysis of bio-derived waste has the potential to provide a lower-carbon, more sustainable way of producing chemicals that are currently made from petrochemicals. But with biowaste feedstocks being so diverse, how can we determine the best use of a particular resource? We talk to Dr Daniel Nowakowski at Aston University's Energy and Bioproducts Research Institute (EBRI) about the challenges faced when thermally processing biowaste, and how collaborating with Shimadzu is helping him and his team to maximise the potential benefits of these feedstocks.

IET JANUARY / FEBRUARY **2024**

Adam Elaradi (Undergraduate Researcher) and Dr Sarah Asplin (Postdoctoral Research Associate) operate the laboratory-scale intermediate pyrolysis system at EBRI. The system can take up to 300 g/h of biomass, and has a maximum processing temperature of 600 °C.

are generally best turned into biochar because of the low loadings of contaminants in the final products, whereas low-quality feedstocks are better suited to bioliquid or biogas, because clean-up by liquid–liquid separation, distillation or downstream processing is a necessary part of producing them".

A relatively new development in processing technology is socalled 'intermediate pyrolysis', whereby the feedstock is placed on a moving bed, she adds. "This allows a high degree of control over the process temperatures and times, making it possible to fine-tune the conditions to optimise the desired products".

Quick initial analysis – Combining the right methods

To properly understand the potential of biomass and other organic waste requires an in-depth study of their chemical composition, says Dr Nowakowski: "You can't simply get away with applying a single analytical method!". In fact, his understanding of this aspect extends beyond his research, because he also has overall responsibility for the wide range of analytical equipment at EBRI.

The advantage of this setup, says Dr Nowakowski, is its speed – "in just an hour, you can complete the preparation and analysis of a test sample, and get an idea of the vapours that could be obtained from the feedstock" he explains. By adding an internal

One of the first steps when understanding the potential of a new material, he explains, is to carry out a thermal degradation study. This involves heating a small portion (about 10 mg) up to about 750 °C in a thermogravimetric analyser and monitoring how the mass changes as the temperature rises, giving the researchers an initial idea of its volatile content and solid residue. At the same time, the gases released can also be sent straight to a detector using a technique known as evolved gas analysis, again to provide a first impression of what the sample contains.

More information on the products obtained from thermal degradation is then obtained by micropyrolysis–GC–MS. This uses what is in essence a miniature version of the fullscale pyrolysis equipment, with a gas chromatograph–mass spectrometer fitted to the outlet, configured to analyse the vapour-phase products. The research team at EBRI have two such 'pyroprobe' systems, with the one they use the most being a multi-shot micropyrolyser from Frontier Laboratories partnered with a Shimadzu GC–MS.

standard, the amounts of given products can be quantified, and a preliminary assessment of product yields obtained, he adds.

The next steps depend on the outcome of those initial investigations, says Dr Nowakowski, but they will often employ elemental analysis, or TOC analysis (using a Shimadzu instrument) to find the 'total organic carbon' present in the sample. Another method useful at this point is calorimetry to find the heating value. "Both these techniques can be useful in discussions with clients who might be wondering whether it's worth simply burning the biowaste to release its energy content", he says.

Detailed analysis of the product mix

When it comes to more detailed analysis, the team turn to a range of chromatographic techniques to cover every possible combination of feedstock type, mixture complexity, and analyte chemistry.

- **• GC–MS/FID,** by splitting the column outflow and sending it two ways, combines the compound-identifying capabilities of MS with the quantitative performance of FID. There are two such systems at EBRI, both Shimadzu – one with a one liquid autosampler for investigating bioliquids, and the other with a headspace-configured 'pyroprobe'.
- **• GC–FID** (gas chromatography–flame ionisation detection) offers simpler and more accurate quantitation than GC–MS, and is used for quick analysis at every stage of biowaste processing. At EBRI, there are a total of six Shimadzu systems that are ready to go with different chromatographic columns plus two systems that also have capability for TCD (thermal conductivity detection). This is used instead of FID when the analytes don't contain a carbon–hydrogen bond, making it

suitable for analysing permanent gases such as hydrogen, carbon monoxide, carbon dioxide, oxygen and nitrogen. One of these GC–FID/TCD systems has a high-temperature injection port for analysing the heaviest hydrocarbons.

• GC–BID is another system at EBRI that's tailored to gas analysis, and uses Shimadzu's BID (barrier ionisation detector), which offers much lower detection limits than TCD for permanent gases.

• **HPLC** (high-performance liquid chromatography) is the standard method for separating gram-quantities of complex liquid mixtures, and there is a Shimadzu system at EBRI used for analysing aqueous phases from the pyrolysis process, especially organic acids.

As is clear from the above list, Shimadzu systems are a big feature of the labs at EBRI, with over 12 instruments in total. "Having the equipment from the same manufacturer means that we can use the same software and the same analytical processes – which is good both for productivity and training", says Dr Nowakowski.

And, says Dr Nowakowski, this equipment and expertise is shared with academic institutions, R&D centres and industrial partners at home and abroad – such as the Biochar CleanTech Accelerator. "Doing this allows us to exchange scientific knowledge, build teams for larger-scale research proposals, and most importantly focus on applied research", he explains: "By carrying out such 'proof-of-concept' studies, we can make big strides towards product certification and commercialisation".

EBRI's impressive array of analytical equipment supports is also available to young researchers, he adds: "We have links with many academic institutions from the UK and Europe – all of whom can apply to use our equipment", he says. "So at any one time, we might be hosting several Masters and Ph.D. students, based here for anything from a few weeks to several months, and getting the benefit from our instrumentation".

Biochar as a soil improver and polymer additive

So what projects is Dr Nowakowski currently involved with?

One project with clear benefits on the ground, he says, is a collaboration with Birmingham City Council. This involves generating a biochar-based soil improver from the large amount of tree and shrub cuttings that the council produces from the parks and gardens that it maintains. Called the 'Urban Biochar and Sustainable Materials Demonstrator' project, the cuttings are taken to the council-run Cofton Nursery in Longbridge (Birmingham, UK), and first screened to remove leaves and other compostable material. The larger more woody material is shredded and pyrolysed on-site in a unit installed in a shipping container, yielding biochar. This is used in various projects locally to improve soil structure and boost plant growth, while also incorporating the carbon in a form that is not readily biodegraded, and so providing a role as a 'carbon sink'.

Nowakowski explains that EBRI was closely involved in optimising this setup: "At the outset of the project in 2019, we investigated the potential for using biochar in wide range of applications. We did quite a bit of large-scale work to find the optimum process parameters to produce larger quantities of biochar", he says.

Another use of biochar that Dr Nowakowski's team is investigating relates to its use as a filler in plastic composites, used (for example) in 3D printing. This is relevant because, as he explains, biochar isn't always 100% carbon: "Depending on the feedstock and process conditions, cold spots can arise in the biochar, which can trap small amounts of low-volatility compounds such as polycyclic aromatic hydrocarbons, which are toxic. These would normally remain bound to the biochar, but if you're using it as a filler in plastic, which then ends up being heated during processing, it could release these chemicals and cause a safety issue".

The analytical framework for pyrolysis of biowaste used at EBRI, showing the main stages (circles) and products (rectangles), and the analytical processes used at each stage.

• Organic chemicals: GC-MS/FID, HPLC

Anitta Xavier (Research Placement Student) injects a biomass sample into a Frontier Lab micropyrolyser for direct transfer to the Shimadzu GCMS-QP2010 SE system.

WWW.**ENVIROTECH-ONLINE**.COM

32 **Environmental Laboratory**

IET JANUARY / FEBRUARY **2024**

To study this effect, he says, the researchers use the 'pyroprobe' setup to assess the thermal stability of small samples of biochar-modified plastic pellets, to be sure that the product is going to be safe for the purpose.

Bioliquids as fuel additives and antifungal agents

Bioliquids are also a fruitful area of study, says Dr Chris Thomas, Knowledge Exchange Associate at EBRI, and another former Ph.D. student of Dr Nowakowski. "A particular interest at the moment is what we call the 'catalytic upgrading' of pyrolysis liquids. This involves taking the crude, water-immiscible pyrolysis oil and treating it with hydrogen under high pressure and temperature with a catalyst. That removes oxygen from the oxygen-containing compounds, giving us an upgraded pyrolysis liquid, which can then be distilled and combined with conventional biorefinery streams for production of biofuels.

For this application, says Dr Thomas, the boiling-point distribution is of particular interest, but distillation doesn't work well on small scales, and optimising the pyrolysis–upgrading process on a large scale would be expensive and timeconsuming. So to predict the boiling-point properties quickly from smaller-scale experiments, he used one of the GC–FID systems at EBRI to obtain a chromatogram of the crude liquid, and then applied Shimadzu's SimDist software tool to interpret the results.

Dr Thomas is keen to emphasise the benefits of this approach: "Directly from what are often very complex chromatograms, we can use this 'simulated distillation' method to help us understand at what temperature the different components in a mixture will boil off at. The resulting distillation curve shows us the proportion of each component that evaporates at different temperatures, which is crucial for checking the quality of biofuels and other bioliquid derived chemicals".

Another application of bioliquids that Dr Nowakowski and colleagues are working on is so-called 'wood vinegar'. This is a light-coloured aqueous liquid dominated by carboxylic acids, including acetic acid, formic acid and propionic acid, as well as some phenolics, and many other organic compounds about which little is known.

Dr Nowakowski explains: "These liquids have received a lot of hype as 'natural' antifungals and antibacterials, but there's little information about exactly what they contain, let alone studies into their environmental and health impacts. So a focus for us at EBRI over the coming year or so will be using our HPLC and GC– MS instruments to understand these liquids in more detail".

This research is vital, he adds. "Wood vinegar definitely has potential for treating mould and bacterial films, but there's a big issue around process and quality control. For example, if the pyrolysis temperatures are too high, the product can contain toxic aromatics, dioxins, and chlorine-containing chemicals. If you're applying the wood vinegar to crops, they'll end up in the environment, which is something you certainly don't want".

Diverse feedstocks, multiple processes

This is just a flavour of the work that's going on at EBRI, says Dr Nowakowski. Other projects involve looking at other sorts of waste, including chicken manure, municipal solid waste, used cooking oil, spent barley grains from beer production, bicycle tyres, offcuts of aluminium–plastic 'thermal covers' and cork, as well as a scoping exercise for carbon sequestration using farmed seaweed.

And it's not all about pyrolysis, either. When the feedstock contains a lot of water that would be too energy-intensive to remove – such as algae – they can use a process called hydrothermal liquefaction. This involves catalytic treatment of a liquid slurry, with the result being a 'biocrude' similar to that obtained from pyrolysis of the solid material.

Biological processes such as anaerobic digestion and fermentation are also researched at EBRI, says Dr Nowakowski. "When a feedstock has very low calorific value but contains some good sugars, then we can get microorganisms to convert it to methane or ethanol, which can then be used as a fuel. The advantage of this approach is that the end product is usually quite refined, making the analysis a lot simpler".

In summary, it seems that EBRI has a solution for every challenge: "Whatever the feedstock type, we can select and tailor the process to extract the maximum value from it!", he says.

Expanding biowaste research – with support from Shimadzu

Through all of his work as a researcher and laboratory manager at EBRI, Shimadzu has been there for him, explains Dr Nowakowski. "Right from 2013 when we were setting up the new laboratories for bioenergy research here at EBRI, I've been able to benefit from their expertise in analytical chemistry", he says.

"They've always been ready to discuss our needs, and in many cases they've made a major contribution to the decision-making process. For example, in one project we needed to analyse permanent gases from our hydrothermal liquefaction unit, and they helped with setting up a GC–BID system – which significantly improved our research outputs compared to the GC–FID system we'd originally been planning".

This collaboration has continued to the present day, he adds, with guidance on getting the most out of the equipment, as well as routine maintenance and timely support. "With so many people booking our equipment, business continuity is very important for us $-$ so it's great that we always have a quick response from Shimadzu!", he says.

And it's clear that there's no let-up in the pace of work at EBRI. Dr Nowakowski says that they're still expanding their operations: "We're increasingly focusing on applied research and novel processing methods, including biological processes for the production of value-added chemicals. We're also becoming very strong in materials science, with emphasis on carbon materials derived from biochar. In addition, very recently EBRI started development of an integrated laboratory-scale centre to explore production of low-carbon hydrogen from biomass".

"There's so much going on – and it's good to know that Shimadzu will be there for us, ready to help with anything I need!", he concludes.

Author Contact Details

Shimadzu UK Limited

• Address: Mill Court, Featherstone Road, Milton Keynes. MK12 5RE

• Tel: +44 (0)1908 552209 • **Web: www.shimadzu.co.uk**

A typical example showing the results of analysing a sample of biochar by GC–MS to assess its purity. In this case, the compounds highlighted would be undesirable in the final product because of their known toxicity; work carried out at EBRI aims to understand the pyrolysis conditions that lead to these contaminants, and minimise them. The analysis used a Frontier Lab EGA/PY-3030 double-shot analytical pyrolyser, close-coupled with a Shimadzu GCMS-QP2010 SE system.

Left: Sachin Solomon (Knowledge Exchange Associate) uses the Shimadzu GCMS-QP2010 SE system with an AOC-20i autoinjector to carry out the GC– MS analysis of pyrolysis bioliquids. Right: Dr Christopher Thomas (Knowledge Exchange Associate) working with EBRI's 'simulated distillation' system, consisting of a Shimadzu Nexis GC-2030 with SimDist software.

A typical example showing the results of analysing a sample of biochar by GC–MS to assess its purity. In this case, the compounds highlighted would be undesirable in the final product because of their known toxicity; work carried out at EBRI aims to understand the pyrolysis conditions that lead to these contaminants, and minimise them. The analysis used a Frontier Lab EGA/PY-3030 double-shot analytical pyrolyser, close-cou-

pled with a Shimadzu GCMS-QP2010 SE system.

RESEARCHER BIOGRAPHY

Dr Daniel Nowakowski is a Lecturer in Chemical Engineering and Applied Chemistry at Aston University (Birmingham, UK), and Laboratory Manager of Aston's Energy and Bioproducts Research Institute (EBRI). Following undergraduate studies in Chemical Technology in Gliwice (Poland), he completed his Ph.D. degree

in Environmental Catalysis at the University of Leeds (UK) in 2007. The same year, he joined Aston University as a post-doctoral research assistant, and since 2012 he has held several senior fellowship posts. He has been an active researcher since 2007, and has been involved in a wide variety of research projects and collaborations relating to the processing, analysis and application of biomass and pyrolysis products.

READ, **SHARE** or **COMMENT** on this article at: **envirotech-online.com/article**

